


How to install OlympusViewer Plugin

Windows


- 1 Download OlympusViewer-win.zip.
- 2 Extract the zip file.
- 3 Execute OlympusViewer-win.exe
- 4 Read END-USER LICNESE AGREEMENT
If you agree to it, click Yes.
If you don't agree it, click No.

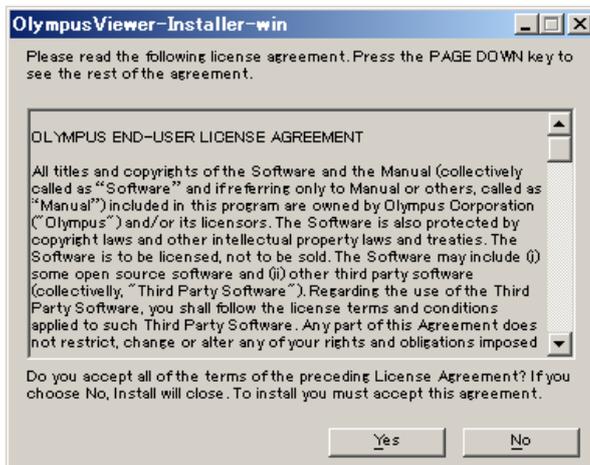

- 5 Specify your extraction path:



- 6 Extract OlympusViewer-package.zip

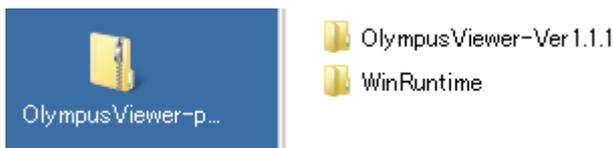

7 Install VS2008 redistribution package:

If you use 32bit os, install vcredist_x86.exe

If you use 64bit os, install vcredist_x64.exe



8 Install OlympusViewer plugin:

Copy OlympusViewer folder to your ImageJ plugin folder.

If ImageJ plugin folder already has OlympusViewer folder, delete the folder before copying.




Mac


- 1 Download OlympusViewer-mac.dmg.
- 2 Double click the dmg file.
- 3 Read END-USER LICNESE AGREEMENT

If you agree to it, click Agree.

If you don't agree it, click Disagree.

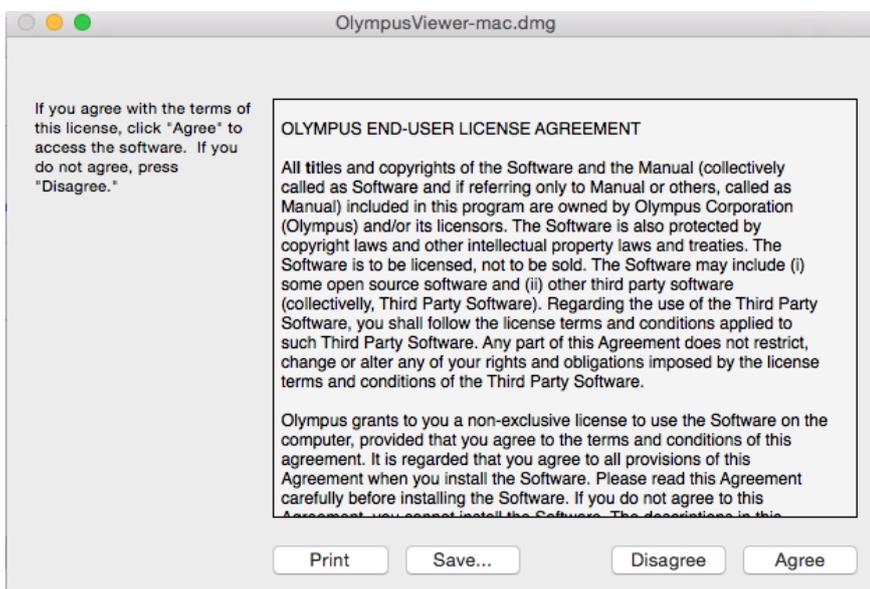

- 4 Install OlympusViewer plugin:
Copy OlympusViewer-Ver2.1.1 folder to your ImageJ plugin folder.
If ImageJ plugin folder already has OlympusViewer folder, delete the folder before copying.

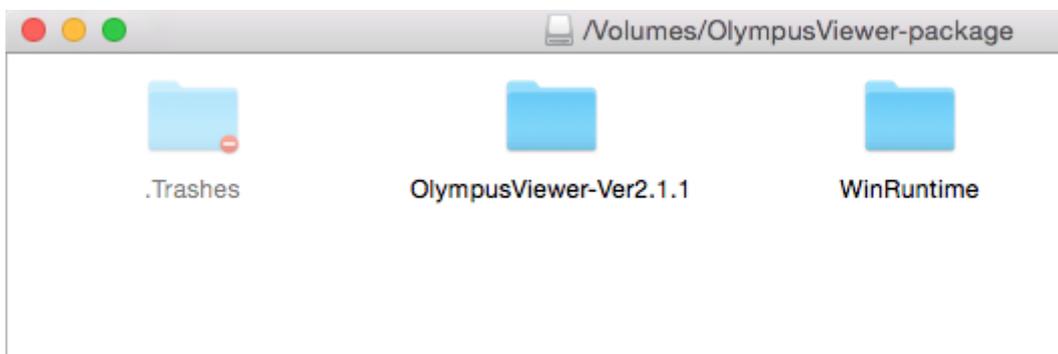

Read END-USER LICNESE AGREEMENT